

Cellovations® Human Hepatocyte Thawing Medium

Cellovations® Human Hepatocyte Thawing Medium

- optimized thawing procedure for all Hepatocytes

Product Description:

Cellovations® Human Hepatocyte Thawing Medium is a complete medium formulated for optimal thawing of all hepatocytes, human and animal, for 2D&3D application. It contains essential and non-essential amino acids, vitamins, organic and inorganic compounds, hormones, trace minerals together with the growth factor containing supplements.

The thawing medium guarantees a high viability of hepatocytes.

The medium is for research use only and not approved for human or animal use or for application in any in vitro diagnostic procedure.

Storage of the medium and supplements

The hepatocyte thawing medium will be delivered in two vessels (Compound I + II). Mix both compounds and use the thawing medium the same day!

Please note that the medium has to be stored at 4°C but never should be frozen.

Preparing and Stability:

Warm up the medium. Please rinse the bottle with 70% ethanol as protection. Remove ethanol excess before opening the cap of the Bottle. The complete medium should be used within four weeks.

Please be careful not to touch the bottle interior threads with your fingers. Please avoid long term exposure to light because some components are light sensitive. Do not warm up the medium over 37°C. All this protects your medium against contamination or degradation of essential components.

For more details how to use the medium see our instructions.

• Catalogue number: PB-MH-053-5900-T-50 | -250 | -500

Components: 50 or 250 or 500 ml Hepatocyte Thawing Medium.

ORDER UNDER sales@pelobiotech.com or FAX + 49 (0)89 517 286 59 88