

Cellovations® Endothelial Cell Growth Media

Cellovations® Endothelial Cell Growth Medium Kit classic-GFP/RFP

- optimized growth conditions for Endothelial Cells;
especially GFP/RFP labeled cells

Product Description

Cellovations® Endothelial Cell Growth Medium Kit classic-GFP/RFP is a complete medium formulated for optimal growth of endothelial cells that were labeled with GFP or RFP. It contains essential and non-essential amino acids, vitamins, organic and inorganic compounds, hormones, growth factors, trace minerals plus FCS together with the supplements.

The medium is for research use only and not approved for human or animal use or for application in any in vitro diagnostic procedure.

Storage

Please note that the basal medium has to be stored at 4°C up to 8°C but never should be frozen.
The supplements has to be stored at -20°C. Protect from light for a longer time period.

Preparing and Stability

Thaw the supplements at 37°C before adding to the basal medium. After thawing please rinse the bottle and the tube with 70% ethanol as protection. Remove ethanol excess before opening the cap of the tube. After adding the supplements to the basal medium slew carefully to get a homogenous medium.

Please be careful not to touch the tube or bottle interior threads with your fingers. Please avoid long term exposure to light because some components are light sensitive. Also do not warm up the medium over 37°C. All this protects your medium against contamination or degradation of essential components.

Product Information

- Catalogue Number: PB-MH-100-2190-GFP/RFP
- Components:
 - 500 ml Basal Medium
 - Supplement Kit
 - Glutamin - FCS
 - bFGF - EGF
 - Hydrocortison - ECGS/H
 - optional: Gentamycin/Amphotericin 50 µg/ml

FOR R&D USE ONLY!

MSDS are available upon request!

ORDER UNDER sales@pelobiotech.com or FAX + 49 (0) 89 517 286 59 88